Development of 3D bioprinted human breast cancer for in vitro screening of therapeutics targeted against cancer progression

Shelby M. King, Vivian Gorgen, Sharon C. Presnell, Deborah G. Nguyen, and Benjamin R. Shepherd Tissue Applications and Systems Engineering Groups

Presentation Number: 1805

Abstract

During their lifetime, 1 in 8 American women will develop breast cancer, and an urgent need exists for targeted, safe therapeutics, particularly for refractive disease. Regulatory agencies are currently seeking improved preclinical oncology models to overcome hurdles to efficient drug development, including bridging the gap between the use of 2D cell lines and 3D animal models. To more accurately model initiation and progression of breast cancer and to more efficiently define targeted therapeutics, the breast microenvironment must be considered. Breast stroma, composed of fibroblasts, endothelial cells, and adipocytes, plays a key role in the process of carcinogenesis and metastasis. These cell types secrete extracellular matrix, growth factors, and hormones that affect how therapeutic agents access and target cancer cells. We have used Organovo's NovoGen BioprintingTM Platform to develop a model of human breast cancer in which breast cancer lesions were surrounded by a physiologically-relevant stromal milieu consisting of adipose (differentiated from mesenchymal stem cells), mammary fibroblasts, and endothelial cells. This system has several advantages over current screening tools, including the ability to simultaneously measure the effects of small molecules on cancer cells as well as different cell types in the breast microenvironment. Histomorphological analyses of bioprinted neotissues demonstrated that they were stable and viable for at least 14 days in culture and characterized by clear compartmentalization of adipose, stromal, and epithelial components. During the bioprinting process, robust microvascular networks were created using endothelial cells as a component of the 3D tissue design. Mammary neotissues were fabricated directly into multi-well plates and used to establish biological response profiles to signal mediators including estradiol, progestin, prolactin, and HGF as well as the standard chemotherapeutic agents cisplatin, paclitaxel, methotrexate, and tamoxifen. The effects of chemotherapeutic drugs on specific cell types within the neotissues was assessed by staining with cell-type specific markers in addition to fluorescent live/dead and cytotoxicity assays. The response of the 3D breast neotissues to growth factors, hormones, and chemotherapeutic agents was compared to the response of 2D breast cancer cell lines. This model can thus be used to screen new anti-cancer targets with better efficiency and accuracy for targeting of cancer cells in the context of the in vivo microenvironment.

The NovoGen MMX[™] Bioprinter



Figure 1. The Novogen MMX[™] Bioprinter deposits multiple types of cell aggregates in spatially-defined patterns. Following bioprinting, the cells secrete extracellular matrix and spontaneously organize into relevant tissue microarchitectures in the absence of exogenous scaffold material.

Organovo, Inc., 6275 Nancy Ridge Dr., Suite 110, San Diego, CA, USA





Figure 2. Left, schematic of bioprinted neotissues. A nodule of human breast cancer cells is surrounded by a stromal compartment composed of endothelial cells, fibroblasts, and adipocytes. Right, bioprinted neotissues immediately after printing. Cancer cells are labeled with blue dye.

Histologic Analysis of Bioprinted Neotissues



Figure 3. Histology of bioprinted neotissues. Tissues were analyzed to determine tissue architecture and relative positions of cell types. In panels C-E, breast cancer cells were labeled with CellTracker Green CMFDA (Invitrogen, Carlsbad, CA).



Figure 4. Breast cancer cell (2D) or neotissue (3D) response to tamoxifen. Cancer cells grown in 2D (A) or bioprinted neotissues (B) were treated with media alone, DMSO, or 10 uM tamoxifen for 48 hours and assessed for viability by ATP luciferase assay. For each graph, * indicates p<0.05 for treatment compared to control.

- cells
- 2. Following bioprinting of the stromal compartment, formation of endothelial networks and differentiation of adipocytes were observed.
- 3. Isolated 2D cancer cells were more susceptible to tamoxifen-induced toxicity than cells incorporated into 3D bioprinted constructs when treated with the same dose of tamoxifen for the same duration.

Any statements contained in this report and presentations that do not describe historical facts may constitute forward-looking statements as that term is defined in the Private Securities Litigation Reform Act of 1995. Any forward-looking statements contained herein are based on current expectations, but are subject to a number of risks and uncertainties. The factors that could cause actual future results to differ materially from current expectations include, but are not limited to, risks and uncertainties relating to the Company's ability to develop, market and sell products based on its technology; the expected benefits and efficacy of the Company's products and technology; the market acceptance for the Company's products, and the risks related to the Company's business, research, product development, regulatory approval, marketing and distribution plans and strategies. These and other factors are identified and described in more detail in the Company's filings with the SEC, including its prospectus supplement filed with the SEC on November 27, 2013, its report on Form 10-Q filed November 8, 2013 and its transition report on Form 10-KT filed with the SEC on May 24, 2013 and our other filings with the Securities and Exchange Commission. You should not place undue reliance on these forward-looking statements, which speak only as of the date of this Current Report. These cautionary statements should be considered with any written or oral forward-looking statements that we may issue in the future. Except as required by applicable law, including the securities laws of the United States, we do not intend to update any of the forward-looking statements to conform these statements to reflect actual results, later events or circumstances or to reflect the occurrence of unanticipated events.



2D vs. 3D Drug Response

Conclusions

. Bioprinted tissues retain compartmentalized structures with interaction between stromal and cancer

Safe Harbor Statement